



Installation of LDM-drivers on Windows 10 operating system:

If drivers for the LDM already appear to be installed, start by uninstalling the drivers from the Device Manager menu.

Disconnect the LDM reader, then proceed as follows:

1. Locate the drivers on the installation media (e.g. DMC User). The drivers can also be downloaded from this location: <http://www.mediafire.com/file/xazrawzeqxf3eh/LDM+Drivers.zip>
2. Install the drivers for your reader model (LDM2207/LDM320/LDM330). Make sure to install the drivers for LDM320/LDM330 using admin rights.
3. Connect the LDM reader. It should now be recognized correctly in the Device Manager menu.